





## How to install Driver and application software SK-Viewer

### How to install SK-Viewer and Setting of the adjusted value

1. Insert the CD-ROM to your computer.
2. Double click SK-Viewer 1.658 in the CD-ROM.
3. Start the setup.exe and click "Yes" button.  
Click "Next" button following the dialogues.
4. Click "Done" button to complete the installation.
5. If there is an icon  in your desktop, it has been done successfully.

\*How to use "SK-Viewer" is in another booklet "SK-Viewer operation manual".

6. Double click icon  SK-Viewer in the desktop.
7. Click "Set" button in the software.
8. "Property" window is displayed, click "Advanced settings" button.
9. Choose "Load/Save" tab.
10. Click "Import" button, choose CD-ROM.
11. Click "SKM-Z200C-PC.reg" or "SKM-Z300C-PC.reg" in the CD-ROM and then Click "Open" button.
12. "SKM-Z200C-PC" or "SKM-Z300C-PC" appears and choose it,  
then the colour become reversed blue.
13. Choose "Read" button. \*Reading time is very short.  
Tick "Default load" and click "Apply" button.
14. Click "OK" button.